


THERMAL SOLUTIONS

Characterization of Polymer Film by TMA Penetration

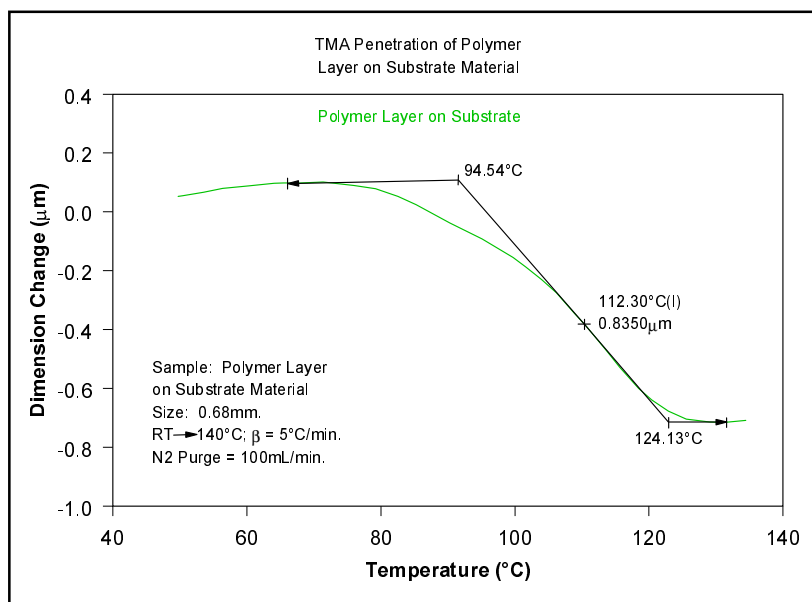

Thermomechanical Analysis (TMA) measures linear or volumetric changes in sample dimension as a function of temperature, time, or force. This data provides valuable information on coefficient of thermal expansion (CTE), penetration, glass transition temperature, softening and flow, and creep/stress relaxation.

This thermal curve shows a TMA penetration experiment on a polymer film bonded to a substrate material. Using a penetration probe, the sample was scanned from room temperature up to 140 $^{\circ}\text{C}$ at 5 $^{\circ}\text{C}/\text{min}$ with a force of 0.1N. A slight expansion of the film is noted before sample penetration begins at 94 $^{\circ}\text{C}$. Total penetration for this sample is measured at only 0.8 μm . The combination of a flat baseline and excellent sensitivity allows the TMA 2940 to measure extremely small changes in sample dimension.

For more information or to place an order, contact:

TA Instruments, Inc., 109 Lukens Drive, New Castle, DE 19720, Telephone: (302)427-4000, Fax: (302)427-4001

TA Instruments S.A.R.L., Paris, France, Telephone: 33-01-30489460, Fax: 33-01-30489451

TA Instruments N.V./S.A., Gent, Belgium, Telephone: 32-9-220-79-89, Fax: 32-9-220-83-21

TA Instruments GmbH, Alzenau, Germany, Telephone: 49-6023-30044, Fax: 49-6023-30823

TA Instruments, Ltd., Leatherhead, England, Telephone: 44-1-372-360363, Fax: 44-1-372-360135

TA Instruments Japan K.K., Tokyo, Japan, Telephone: 813-3450-0981, Fax: 813-3450-1322

Internet: <http://www.tainst.com>

e-mail: info@tainst.com