



TA Instruments

What's New in TRIOS Software

TRIOS Software Overview

TRIOS is TA Instruments' state-of-the-art software package that uses cutting-edge technology for instrument control, data collection, and data analysis for thermal analysis and rheology instruments. The intuitive user interface allows you to simply and effectively program experiments and move easily between processing experiments and viewing and analyzing data. TRIOS software delivers a whole new experiment experience.

- Easy organization and data file management
 - A unique file-naming system allows for effortless organization of data files
 - The File Manager offers simple data file location
- Compatibility with the latest Windows Operating System platforms
- Maximum flexibility
 - Instrument control and data analysis via any computer
 - Configurable for multiple monitors
 - Ability to control multiple instruments at once
- Seamless integration between instrument control and data analysis
- Easy data export in a variety of output formats, including XML, Word, Excel, and PDF
- Effortless graph formatting
- Customization of the display

The TRIOS software is supported by a full range of services, including onsite training, customer service that is only a phone call away, and easy-to-use, easy-to-understand online help. All of these items reflect TA Instruments' commitment to providing thermal analysis and rheology products and related services that deliver maximum value for your investment.

TA Instruments

What's New in TRIOS Software

What's New in TRIOS Software V4.4

TRIOS software is now better than ever with key enhancements that will make your TRIOS experience exceptional. The next generation of instrument control and data collection and analysis, TRIOS V4.4 is more efficient and intuitive, allowing you to work faster and easier.

GENERAL TRIOS ENHANCEMENTS

TRIOS Guardian

Continued development on Guardian technology includes the following updates:

- Added the ability to customize user groups
- Support for file versioning
- Added the option Force Password On Startup
- Added the option Force User Comments on All Results Log Entries
- Added the option to print Guardian settings
- Added Import/Export for easier configuration of Guardian on multiple PCs
- Added the ability to assign Windows groups to Guardian user levels


Localization

Languages have been updated to support Discovery Series DSC 25XX, TGA 55XX, SDT 650, and DMA 850.

RHEOLOGY

Discovery DMA 850

- Miscellaneous bug fixes

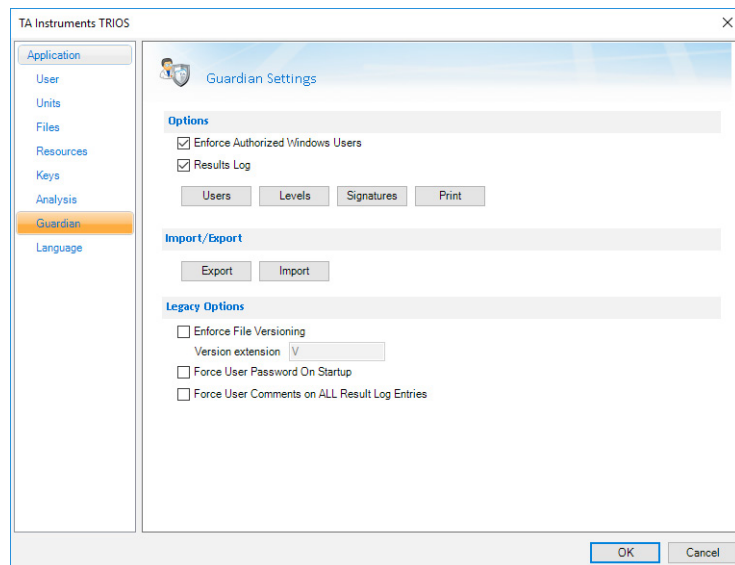

TRIOS Guardian Options

TA Instruments

What's New in TRIOS Software



Additional Information

INSTALLING TRIOS SOFTWARE

For instructions on installing TRIOS software, refer to the [Installing TRIOS Software](#) instructions.

MORE RESOURCES

A number of additional resources are available to you. For assistance with the TRIOS software, first consult the Online Help.

For immediate assistance contact the TA Instruments Hotline at +1 302-427-4000 from 8:00 am to 4:30 pm EST.

For email support, please send your question to one of the following:

thermalsupport@tainstruments.com

rheologysupport@tainstruments.com

microcalorimetersupport@tainstruments.com

PREVIOUS WHAT'S NEW DOCUMENTS

For Previous What's New in TRIOS Software documents, [click here](#).

TA INSTRUMENTS OFFICES

For information on our latest products, contact information, and more, see our web site at: <http://www.tainstruments.com>

TA Instruments — Waters LLC

Corporate Headquarters
159 Lukens Drive
New Castle, DE 19720
USA

Telephone: 302-427-4000

Fax: 302-427-4001

Email: info@tainstruments.com